

Machine learning classification of IMS spectra by **MaSaTECH Chemometrics**

Step1: Prepare classification data

Prepare classification data

Averaged spectrum 2D 2D Selections

Number of classes:

Arabica folders N 19 spectra folders selected.

Robusta folders N 6 spectra folders selected.

Unknown spectra folders 41 spectra folders selected.

Output folder C:/Users/IMS/Desktop/SingleIMS/2DMaps/coffee/Output

Create classification files

Choose if you want to classify : Averaged spectrum along retention time, Complete 2 Dimensional Map or Selected areas from 2D Map

Select number of classes you want to use in your model

Choose recorded spectra of each class. This will create **Training Data**

Choose recorded spectra you **want to classify** (unknown class)

Choose Output folder where to save **prepared data**

Create classification files

Machine learning classification of IMS spectra by **MaSaTECH Chemometrics**

Step2: Principal Component Analysis

Principal component analysis

Averaged spectrum 2D 2D Selections

Num. of components: 20

Decomposition method: Quic

Classification data folder: C:/Users/IMS/Desktop/SingleIMS/2DMaps/coffee/Output

Create PCA files

Choose if you want to use **PCA** for : Averaged spectrum along retention time, Complete 2D Map or Selected areas from 2D Map

Write number of **Principal Components** you want to use for classification




Choose decomposition method for PC generation

Choose folder where you saved your prepared data

Press to create PCA files

Machine learning classification of IMS spectra by MaSaTECH Chemometrics

Step3: Classification

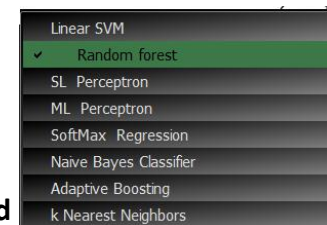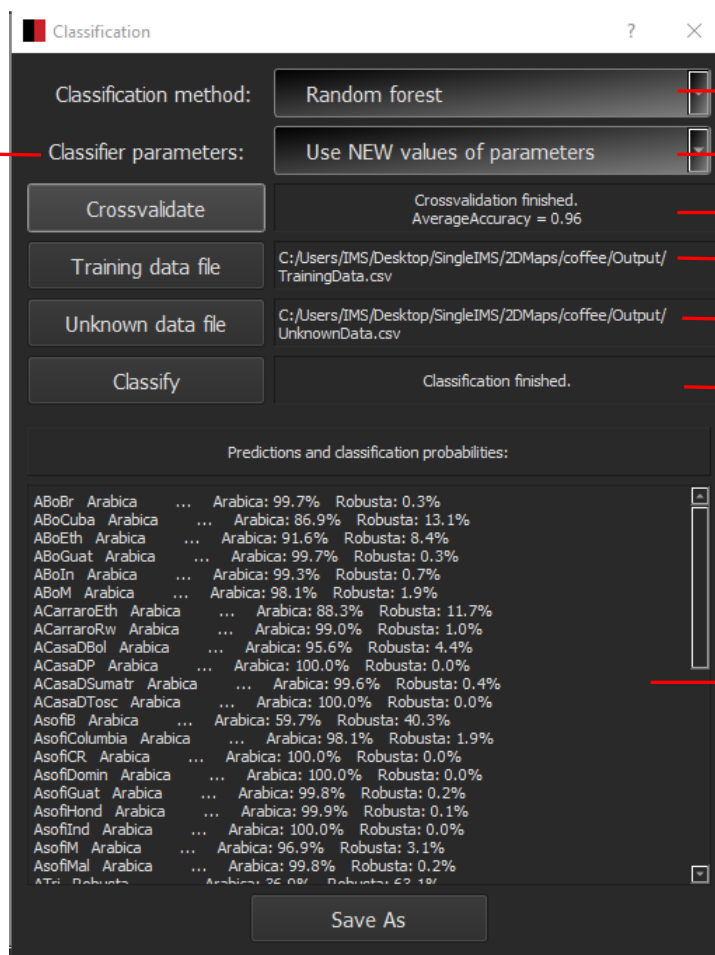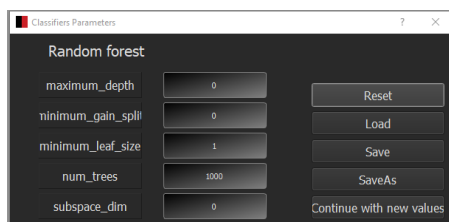

Choose classification method

User have possibility to **modify parameters** of classifier

Calculate **average accuracy** of your model

Select **Training data file** created in step 1

Select **Unknown data file** created in step 1

Press for **classification**

Results gives class information and percentage affiliation to each group.